

Fact Sheet

Instructions for Properly Removing Contaminated Gloves



Figure 1: Grip one glove near your wrist.



Figure 2: As you pull the glove over your fingers, the glove will be pulled inside-out.

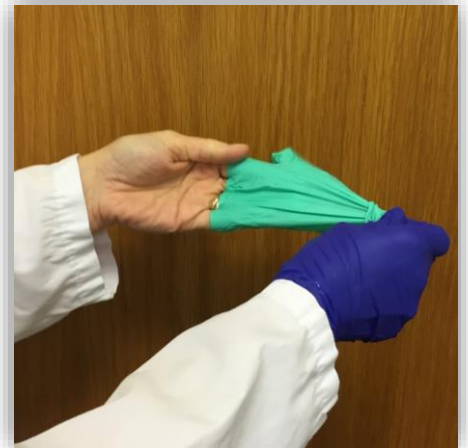


Figure 3: Pull the glove all the way off and grip it in your gloved hand. Be sure to not snap the glove when removing.



Figure 4: Slide your finger under your second glove, still gripping the first glove.

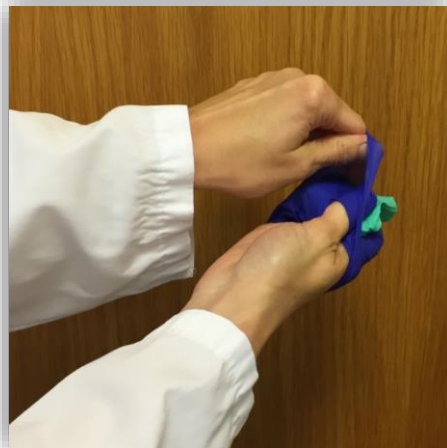


Figure 5: Pull the second glove upwards and off the tips of your fingers, concealing the first glove inside.



Figure 6: Gripping the inside of the removed second glove, the first glove will be contained inside. Dispose of your used gloves properly.